



How to prepare slides for automatic scanning with Panoramic scanner (former Mirax)

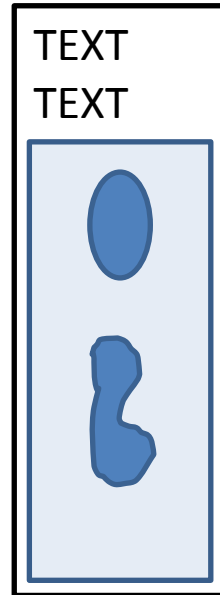

Good

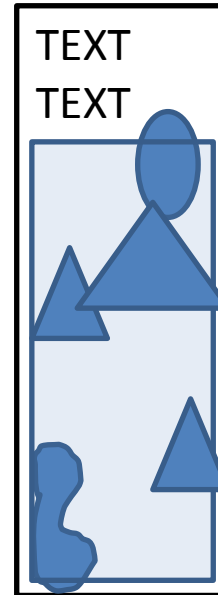

Bad

- Place the sections in the middle of the slide – avoid edges
 - Use coverslips of maximum 50x22mm. Use as little of embedding medium as possible.
- For transmitted light (histo), samples have to be mounted permanently - **slides MUST dry 48h at ca 50°C, or 120h at RT!!!**
- For fluorescence please use curing medium – Dako Fluorescence Mounting Medium **S3023** (almost no air bubbles formed during curing), ProlongGold Diamond, VectaShield Hardset.